



Document No.: BMI_epifluorProtocol_V4[4]

Title: NEON Microbial Productivity from Aquatic Samples

Effective Date: 11/12/2019

Sign-off By

Signatory Role

Sign-off Date

Amy Hill

Document Approver

10/29/2019 6:20:32PM GMT

Janae L Csavina

Document Approver

10/29/2019 9:30:43PM GMT

Lee F Stanish

Document Approver

11/12/2019 11:30:01PM GMT

Jane E Smith

Document Coordinator

11/19/2019 1:46:04PM GMT

NEON Microbial Productivity from Aquatic Samples

Prepared By:
Battelle Memorial Institute
505 King Avenue
Columbus, Ohio 43201

BATTELLE

I. Version History

Previously the percent difference in enumeration (PDE) consisted of monthly analysis of a quality control image by technical staff reporting data. The current version of the SOP describes a PDE procedure where each month 10% of samples (one image per selected sample) will be counted by two technical staff using automated assisted image analysis to verify the Image Analysis process.

II. Objective

The purpose of this protocol is to enumerate microbial cells in water from various domains of NEON within the United States using propidium iodide (PI) staining, black filter filtration, and epifluorescence microscopy. Data from this protocol will be used to determine microbial productivity from sites across the United States.

This protocol is an on-slide staining technique adapted from Boulos et al.¹; modifications include change of fluorophore and concentration, mounting media and volume of sample filtered.

III. Materials and Equipment

Material	Manufacturer	Catalog #
Propidium Iodide (PI)	Molecular Probes	P21493
25mm Black polycarbonate membranes	Sterlitech	1270059
Prolong Diamond	Invitrogen	P36961
Immersion oil		
Coverslips	Premier Brand	SC-92222-1
Microscope slides		
Epifluorescence microscope	Zeiss Axio Imager A.2	N/A

¹ Boulos et al. J. Microbial Methods 37(1):77-86 (1999)

- Milli-Q water (particle-free or 0.22 μm filtered water)
- 30 μM PI working stock
- Fume Cabinet
- Vacuum manifold and pump
- Oven (to dry glassware)
- 25 mm glass vacuum filter holders (including fritted glass funnel and base, aluminum clamps and stopper)
 - All glassware must be particle free
 - Wash glassware using Alconox, rinse in copious amounts of Milli-Q water and dry in oven prior to use

IV. Procedure

A. Sample Filtration and Staining

1. Allow working staining solution (30 μM PI) and Prolong Diamond Antifade Mountant media to reach room temperature. **The stain and ProLong are light sensitive** and should be warmed in the dark (or covered completely with foil).

2. Label microscope slides as follows: **name of sample, date of filtration, analyst initials**

3. Soak black filters in MilliQ water for 1-5 minutes prior to use. Take note of which orientation the filters are in when placed for soaking. The shiny vs non-shiny side of the filter will not be able to be identified after soaking.

4. Place filters shiny side up on the filter manifold. Add 3-5mL of MilliQ water to each filter to both ensure filters stay moist, as well as to check for any leaks in the filter apparatus.

5. *Critical Step:* Filter sample

- a. Add appropriate volume of each sample (or dilution). This should be recorded and is needed to calculate microbial productivity.
- b. Turn on vacuum and slowly filter sample at low (<10 psi) vacuum until all liquid has been pulled through the filter. **Filter to near-dryness, but do not dry the filter – this will result in poor images.**

6. Stain Sample

Note: Steps 6b-6g must occur in the dark. Make sure to work in a dark fume hood or dark area on the bench.

- a. Gently remove filter from the filter manifold with forceps and lay the filter face up on a microscope slide.
- b. Place slide in a dark humid box (sealable box lined with damp paper towels/kimwipes).

- c. Gently pipette the PI working solution over the surface of the filter (filter area should be completely covered with staining solution). Place cover on humid box. Record volume of PI solution used.
- d. Allow stain to sit on the filter for **10 minutes +/- 1 minutes** in the dark.
- e. Stop staining by placing the slide on a dry kimwipe that has been folded over itself a few times. The slide should be placed on the kimwipe at a slight angle, just enough for the solution to soak gently into the kimwipe.
- f. Immediately place 45 µL of ProLong Diamond Antifade Mountant onto the center of the filter and slowly place a coverslip over the mountant media to avoid trapping air bubbles. Do not slide the coverslip relative to the slide or press the coverslip to remove air bubbles.
- g. Allow ProLong to cure for 24 hr at room temperature (RT) in the dark. Slides should be stored at -20 °C in slide box for long term storage after 24 hours.

6. All sample names, dates of filtration, person filtering, filtration volumes, and notes concerning filtration or staining should be recorded

B. Capture Image from Stained Slides

1. Observe slides within 5-7 days.
2. Take images on microscope using a 40X or 100X (preferred) objective lens and save in appropriate data file folder. Image fields should be randomly selected from across the filter. Take at least 15 images per slide. Name images such that they can be tracked back to the specific slide they are acquired from.

Note:

- Counting will be from recorded images, not through the oculars. Analysis of samples is based on the images, so it is imperative to take images in which the number of in-focus cells is maximized.
- When taking images, use the fine focus adjustment to maintain focus – but make certain that all images are taken from the same focal plane.
- Images should be of organisms on the filter and not those in between the filter and the coverslip.
- Avoid collecting image fields off the edge of the filter and around any air bubbles. If a sample has many blank areas or air bubbles, it may indicate a filtration problem and should be noted

C. Assisted Automated Image Analysis using ImageJ

Note: Bacteria densities will range, but generally should be between 5-250 cells per field.

1. Open images using ImageJ software
2. Examine the focus of each image per set

- Select images that are in focus and can be easily adjusted using contrast and brightness.
3. In ImageJ, adjust the parameters of each image for consistent background by selecting “Image” → “Adjust” → “Brightness and Contrast”
 4. Select a contrast range and annotate it on the analysis paperwork
 - a. Have multiple images from the same set opened simultaneously
 - b. Check the setting and ensure that the background staining is eliminated uniformly from each image to be counted.
Note: *the contrast adjustment will vary between different samples and magnifications but SHOULD be the same on all images taken for each slide and magnification.* Be sure to adjust each set before beginning analysis
 5. In toolbar, select “Image” → “Adjust” → “Color Threshold”
 6. Adjust the Brightness setting and annotate the value on the analysis paperwork (Adjusting the brightness setting maximizes the number of cells that ImageJ will recognize).
 - a. Observe each image and adjust each set consistently.
 - b. Do not adjust the Hue setting. This will only change intensity of the color and not the threshold.
 7. Once all appropriate adjustments are complete, select “Analyze” → “Analyze Particles”
 - a. Type in the range 20-2000 as the particle size range to be counted. This is a blanket approach to limit the number of objects identified as microbial cells to those between 20 and 2000 pixels with no relationship to magnification.
 - b. Select “Show Overlay Masks,” “Summary,” and “Display Results”
 - c. The number of counted particles will appear in the “Summary” tab and individual areas are tabulated in “Display Results”

Required if an image has a scale bar: When looking at the image overlay, be sure check the scale bar and omit it from the analysis. This is necessary as the image scalebar is counted as an object when performing the analysis.

8. Record all particles that the software counts and manually count any particles that the software has missed. Record ImageJ counts in the ‘Computer’ column of the analysis spreadsheet.
 - a. In the Threshold Color window, click the Original button to remove the overlays that were not counted by ImageJ.
 - b. Move overlay masks using the computer mouse to determine if there are multiple cells under any of the masks. Count all additional cells not counted by the software.
 - c. Record manually counted cells in the ‘Additional’ column of the analysis spreadsheet.

9. Repeat steps 3 through 8 for all images for each slide across all domains analyzed.

E. Determine Microbial Productivity

Calculate Microbial Cell counts using data captured in step 4 and the following formula:

$$Cf = \frac{(\text{Average number of cells/field}) \times R}{Fs}$$

Where:

Cf = average number cells/mL

R = active area of filter/area of field counted

Fs = volume of initial sample filtered

- For example, if 1 mL of sample is filtered, Fs=1, if 200 µL of sample is filtered, Fs = 0.2
- Fs must consider any dilution made during sample processing for an accurate number to be calculated

Filter Size	Active Filter Area (cm ²)	Area of Image Counted (cm ²)	Objective Used	Conversion factor (R) ^{***}
25 mm	2.1	4.86E-05*	100X	43,209.88
	2.1	8.60E-05**	40X	24,420.51

* Images are 80.56 µm x 60.36 µm using the 100X objective

** Images are 107.13 µm x 80.27 µm using the 40X objective

*** R factor calculated for a specific Zeiss Axiovision A2 Epifluorescence Microscope and color AxioCam MRc5 camera and calibrated Axiovision software. R factor must be calculated for each microscope and camera used for image capture and analysis.

F. Quality Review

To verify the image analysis procedure, 10% of samples each month will be analyzed by two technical staff and the percent difference in enumeration (PDE) will be calculated for each sample. The PDE is the percent difference of the cell counts from each analyst.

PDE determination:

1. 10% of the samples each month will be chosen for quality review (one image per sample selected for the PDE check(s)).

2. Two technical staff each will analyze the images chosen for quality review using the previously described analysis method (Section C, steps 3 through 9). Each technician will be assigned Analyst 1 and 2 prior to doing analysis.

3. The PDEs for each sample will be calculated using data captured in previous step via the following formula:

$$PDE = \left(\frac{|Count1 - Count2|}{Count1 + Count2} \right) * 100$$

Where:

Count 1 = Analyst 1 cell counts

Count 2 = Analyst 2 cell counts

4. If the PDE is <5% (0.05), image counts by the staff member are accepted.

5. If PDE is outside of the acceptable range, the two staff originally performing the PDE check will repeat the count(s). If the PDE continues to be outside of the acceptable range, a third staff member will perform the count(s) and be compared to Analyst 1. Any issues concerning the PDE will be discussed with the Task Lead (TL) before any work continues. The TL will determine if the issues with the PDE check warrant a recount of samples analyzed in the same set as the problem PDE.